

USB DIGITAL CAMERA MICROSCOPE

Users Manual

Installing H-CAM Device Operates Manual

2009/07/27 Update Ver.1.0

Installing H-CAM Device Operates Manual

Contents

1. Please insert CD in CD player	3
2. Install H-CAM application program	4
2.1 Start setup program	4
2.2 Welcome to use H-CAM	4
2.4 Now installing program	5
2.5 Complete install	6
2.6 After installation complete	6
3. Install H-CAM USB digital video driver	7
3.1 Start to install driver program	7
3.2 Welcome to use USB video device setup wizard	8
3.3 Now installing	8
3.4 install complete	9
3.5 After installation complete	9
4. Uninstall H-CAM AP	10
4.1 Choose H-CAM AP's Uninstall driver as Fig. 4-1	10
4.2 Make sure if you proceed it as Fig. 4-2	10
4.3 In process of Uninstall Fig. 4-3	10
4.4 Finish Uninstall	11
5. Uninstall H-CAM USB digital video driver	12
5.1 Choose H-CAM USB digital video driver to uninstall as Fig.5-1	12
5.2 Fig. 5-2 in preparation	12
5.3 If you want to completely remove the driver, as Fig. 5-3	12
5.4 In process of uninstall	13
5.5 Finish uninstall H-CAM USB digital video driver	13

UM-CAM Installing

There are 3 ways to complete H-CAM Installment. Please follow the 3 steps to complete the installation.

- **Step 1:1. Please insert CD.**
- **Step 2:2. Install H-CAM application program.**
- **Step 3:3. Install H-CAM driver.**

There are 2 ways to complete H-CAM Uninstall. Please follow the 2 steps to complete the Uninstall.

- **Step 1: 4. Uninstall H-CAM AP.**
- **Step 2: 5. Uninstall H-CAM USB digital video driver.**

1. Please insert CD in CD player

Please insert the product CD in the player and the computer will read automatically. It will pop up a window shown below:



Fig.1-1 setup window

When the screen can't pop up this window, please go to CD-ROM to select Autorun.exe execute file, which the same will pop up as Fig.1-1 setup window. Autorun.exe is shown below:



Fig.1-2 execute icon

2. Install H-CAM application program

Please select the image 1-1 and setup application program button to install H-CAM program.

2.1 Start setup program

Click setup application program, after it will pop up (please select a language) window on Fig. 2-1, the initial language will follow up the pc OS system. If you want to cancel setup please click“Cancel”then click“OK”.



Fig. 2-1 setup language window

2.2 Welcome to use H-CAM

After choosing the language, go to welcome to use H-CAM setup wizard window fig.2-2, then please click“Next”then continue. If want to cancel please click “Cancel” button..

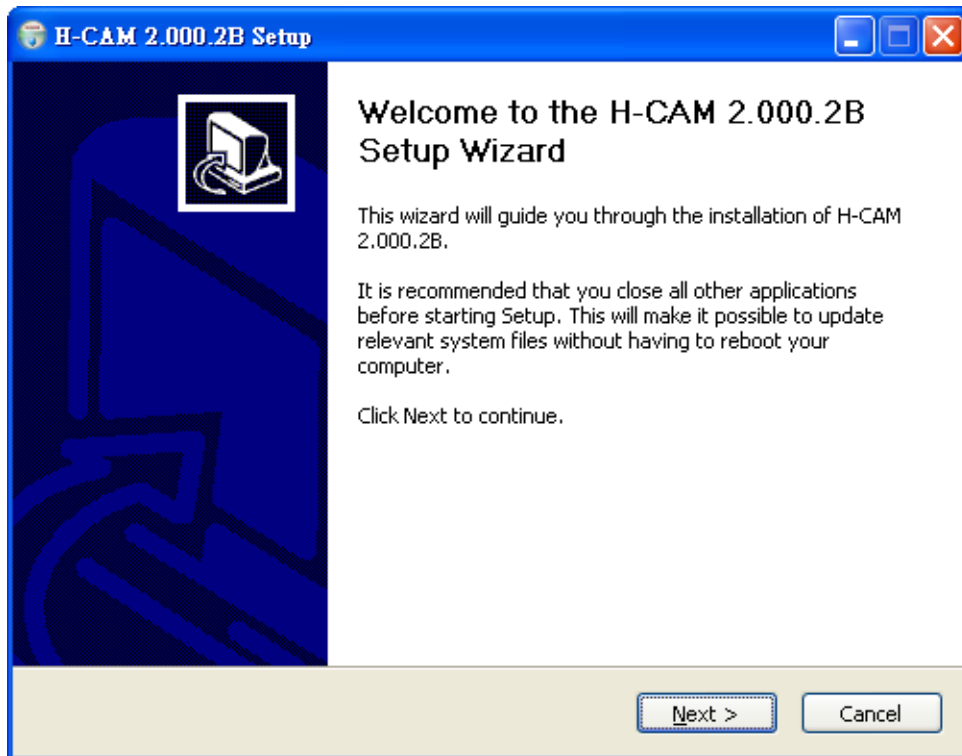


Fig. 2-2 welcome to use H-CAM setup wizard

2.3 Choose installation route

Next step enters choose of installation route window as Fig.2-3. If want to change file route click“Browse”to change setup path, then please click “install” then start installing.

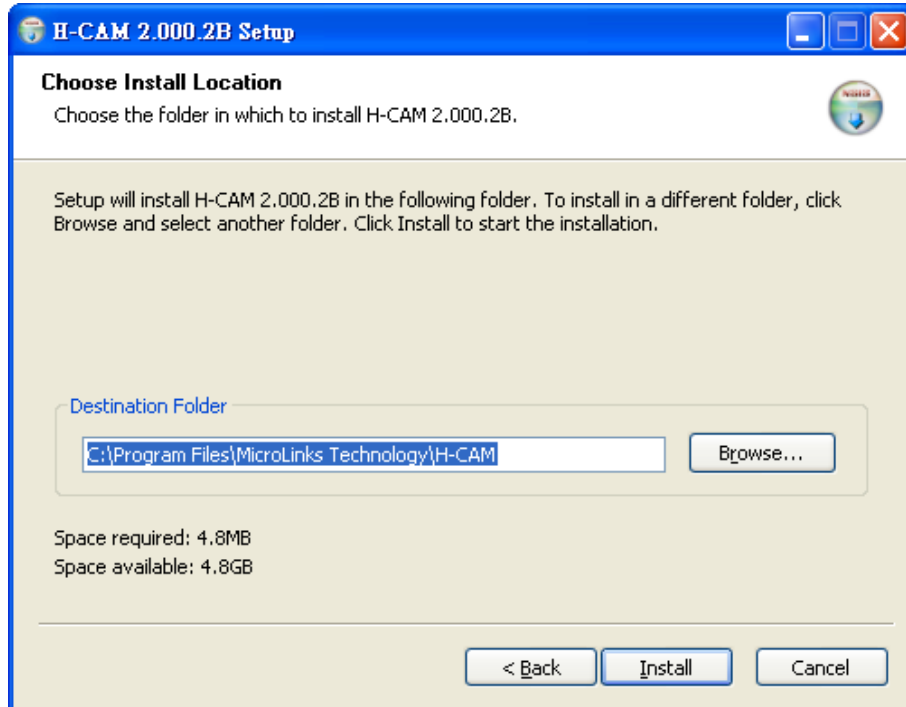


Fig. 2-3 choose installation route

2.4 Now installing program

Continue install wizard will start installing program as directory fig.2-3.

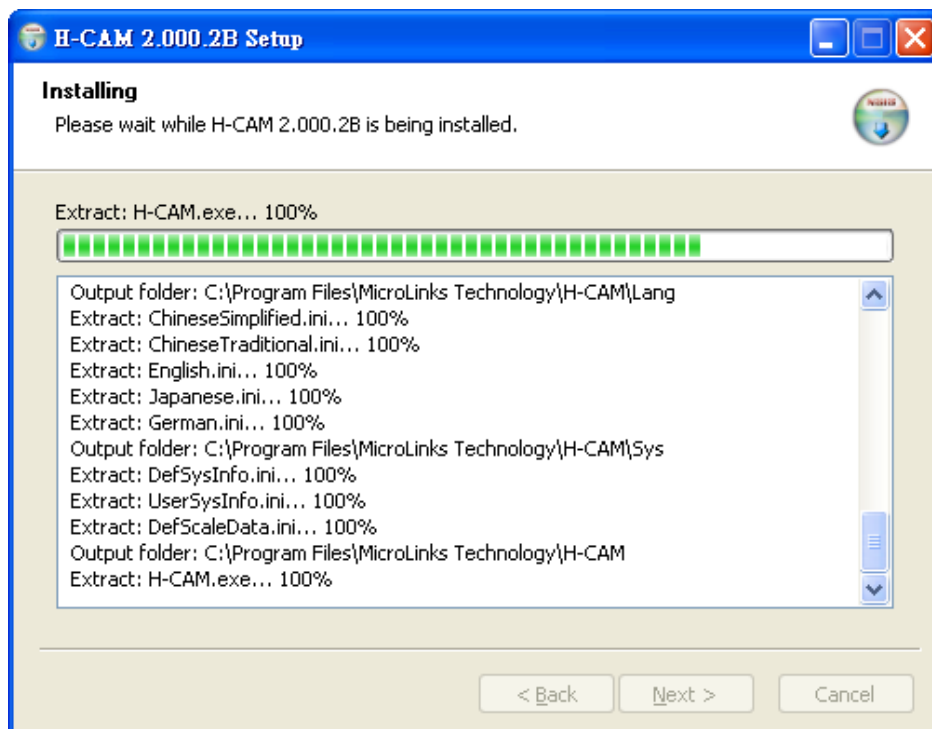


Fig.2-4 now installing

2.5 Complete install

When the H-CAM program complete installation (Fig.2-5). Please click“Finish”.

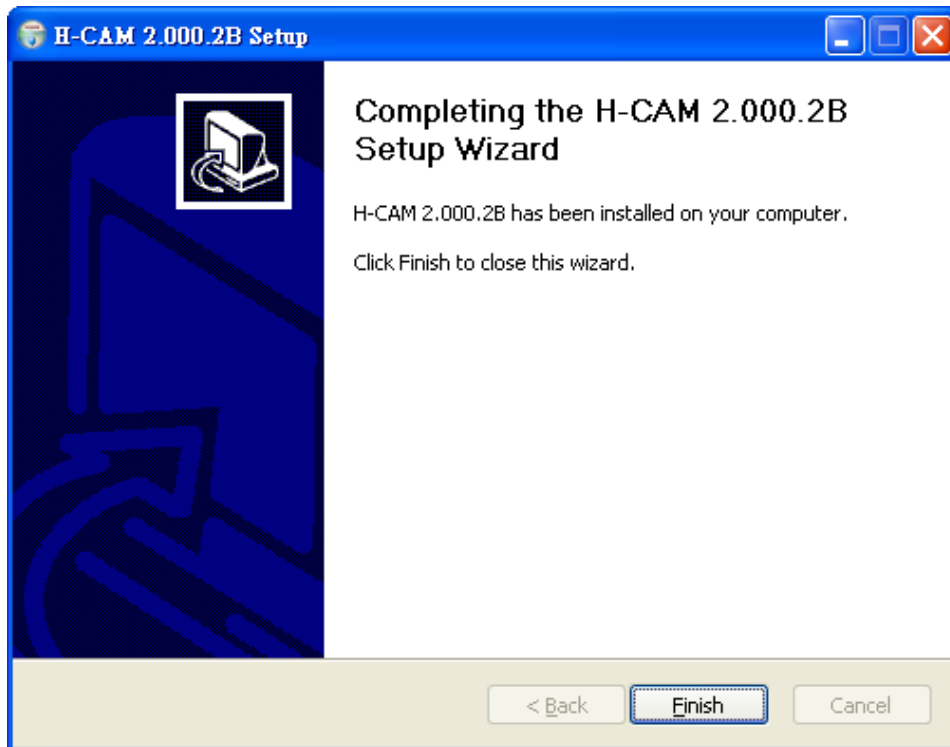


Fig.2-5 H-CAM wizard install complete

2.6 After installation complete

Desktop will show H-CAM short cut as Fig.2-6.

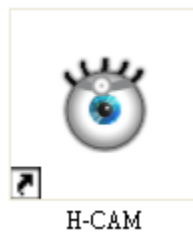


Fig.2-6 H-CAM short cut

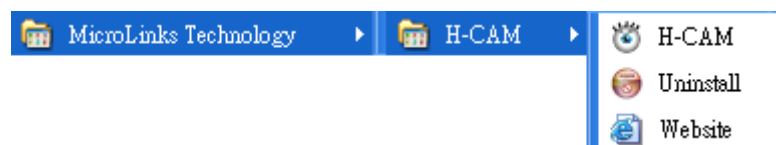


Fig. 2-7 H-CAM at position all programs

3. Install H-CAM USB digital video driver

The driver is optional one, if you install it, you can have more additional functions added to the H-CAM AP.

Attention: Before install driver, please connect H-CAM USB Microscope Camera to the computer. If not, you can't install the program.

3.1 Start to install driver program

Please select Fig.1-1 to start Installing H-CAM drive. Firstly please make sure to connect H-CAM to a pc. Otherwise, it will pop up Fig 3-1.



Fig.3-1 not found H-CAM step up

After click setup H-CAM driver button, it will pop up Fig.3-2, now be ready to install the window. If cancel installation, please click“cancel”button.

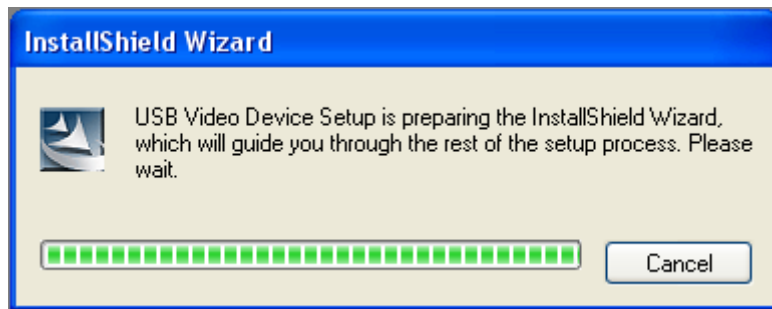


Fig.3-2 Getting ready to install

3.2 Welcome to use USB video device setup wizard

Ready to install, please follow up the steps.

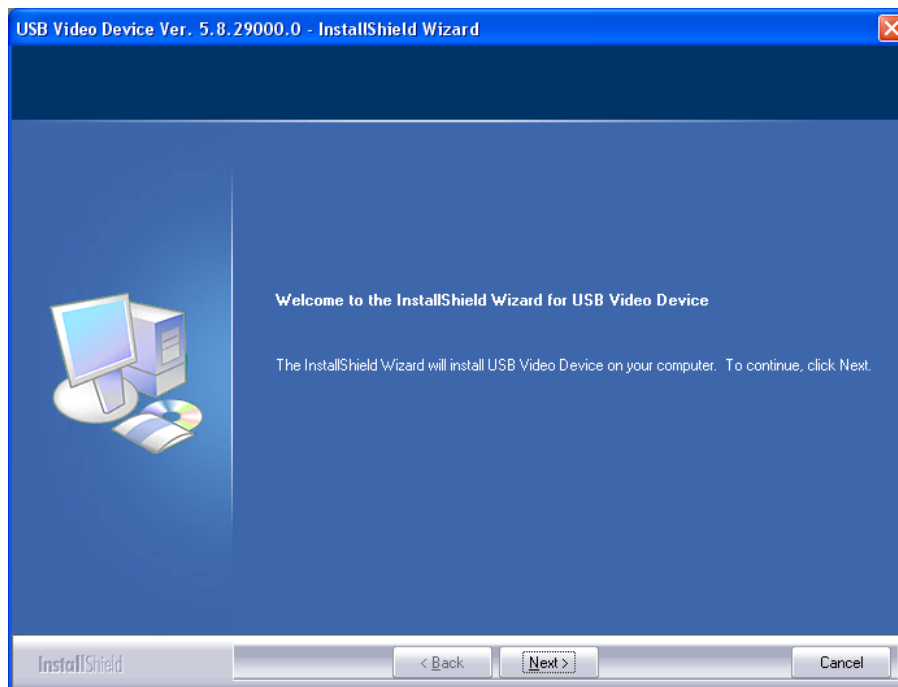


Fig.3-3 welcome to use USB video device Install wizard.

3.3 Now installing

If you want to stop installation, please click“Cancel”.

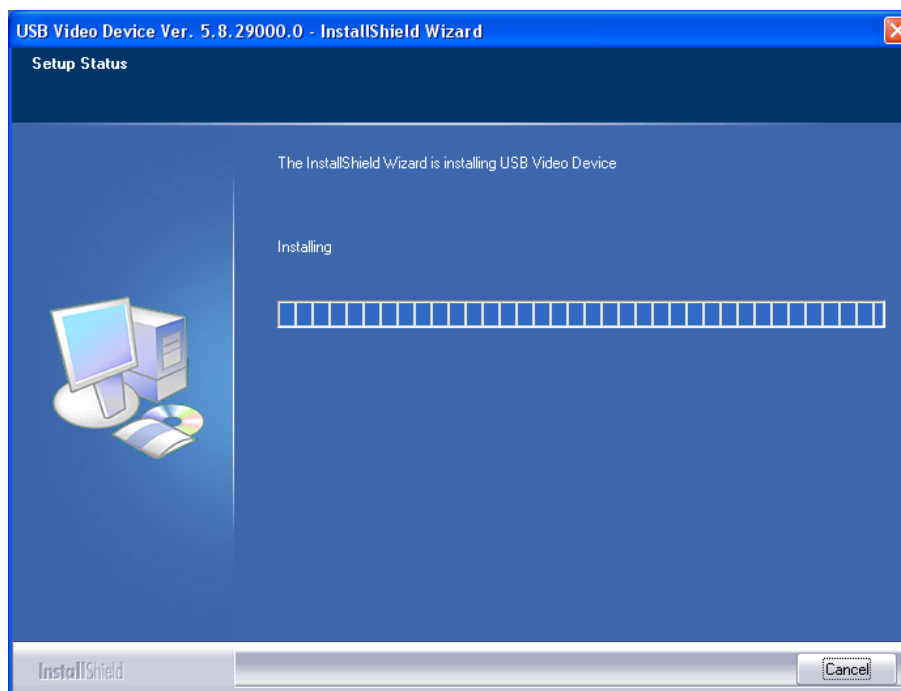


Fig.3-4 Now installing USB video device

3.4 install complete

Complete USB Video Device setup (Fig. 3-5), click “finish” and then driver program is completed.

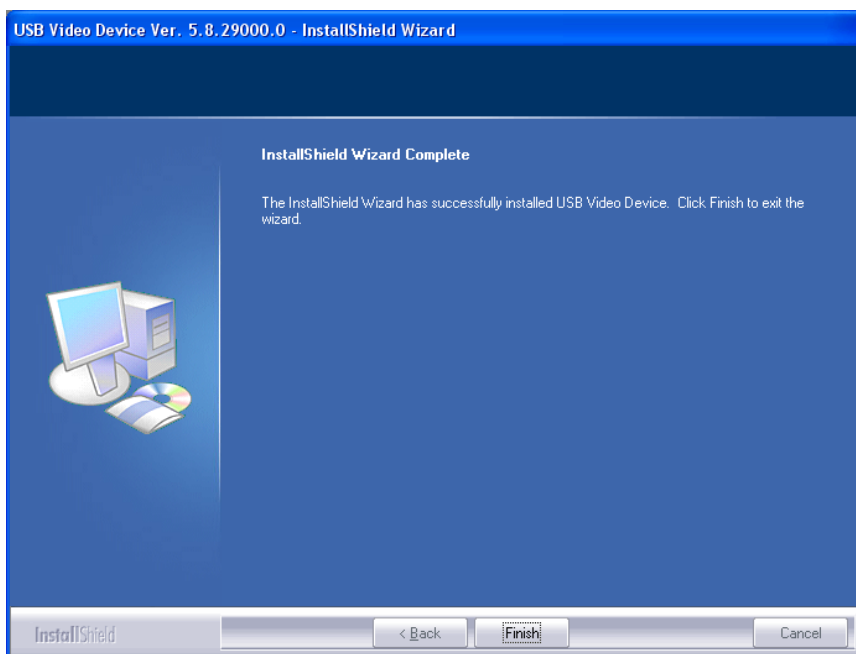


Fig. 3-5 USB Video Device installation complete

3.5 After installation complete

After installation is completed, the window on the right corner will show a video icon Fig.3-6. Start -> all programs will also have USB Video Device file shown Fig.3-7.



圖 3-6 USB Video Device 圖示

H-CAM can be executed from ICON AMCap as below Fig. 3-6. However, AMCap is a simple tool for preview only and can't support image process as the H-CAM AP.



Fig.3-7 USB Video Device is at all programs position.

4. Uninstall H-CAM AP

4.1 Choose H-CAM AP's Uninstall driver as Fig. 4-1

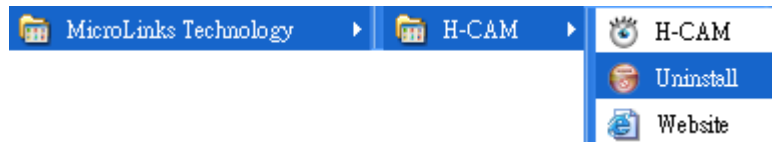


Fig. 4-1 H-CAM AP Uninstall

4.2 Make sure if you proceed it as Fig. 4-2

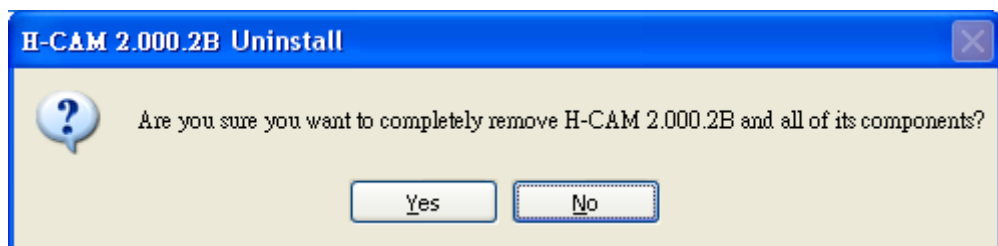


Fig. 4-2

4.3 In process of Uninstall Fig. 4-3

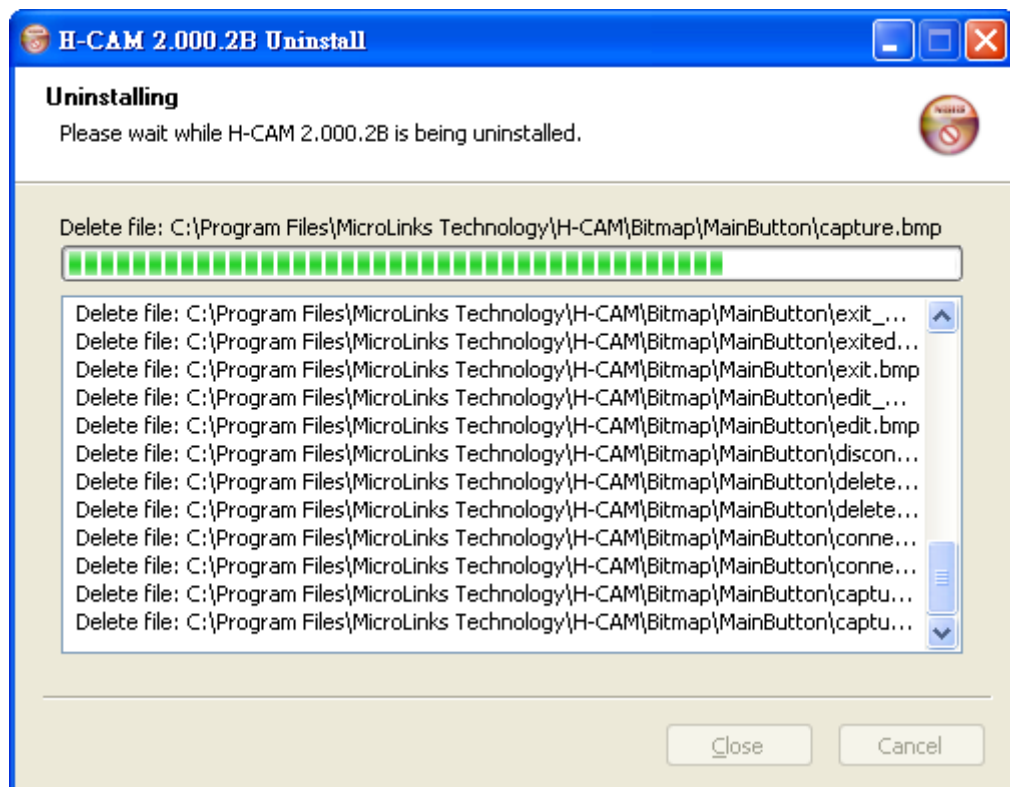


Fig.4-3 in process

4.4 Finish Uninstall



Fig. 4-4 Finish Uninstall

5. Uninstall H-CAM USB digital video driver

5.1 Choose H-CAM USB digital video driver to uninstall as Fig.5-1



Fig. 5-1 U Uninstall H-CAM USB digital video driver

5.2 Fig. 5-2 in preparation

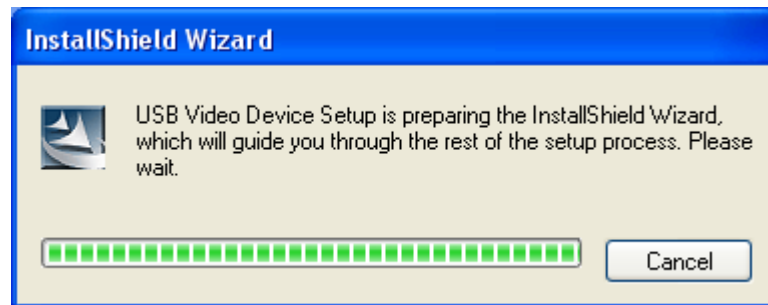


Fig. 5-2 preparation

5.3 If you want to completely remove the driver, as Fig. 5-3.

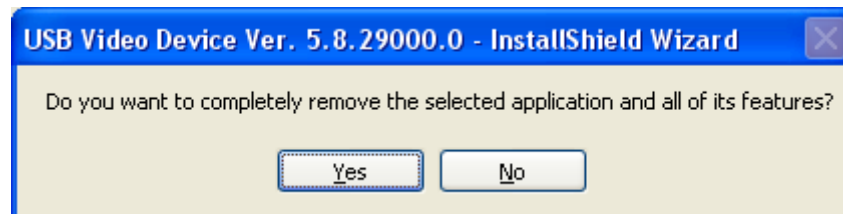


Fig. 5-3

5.4 In process of uninstall

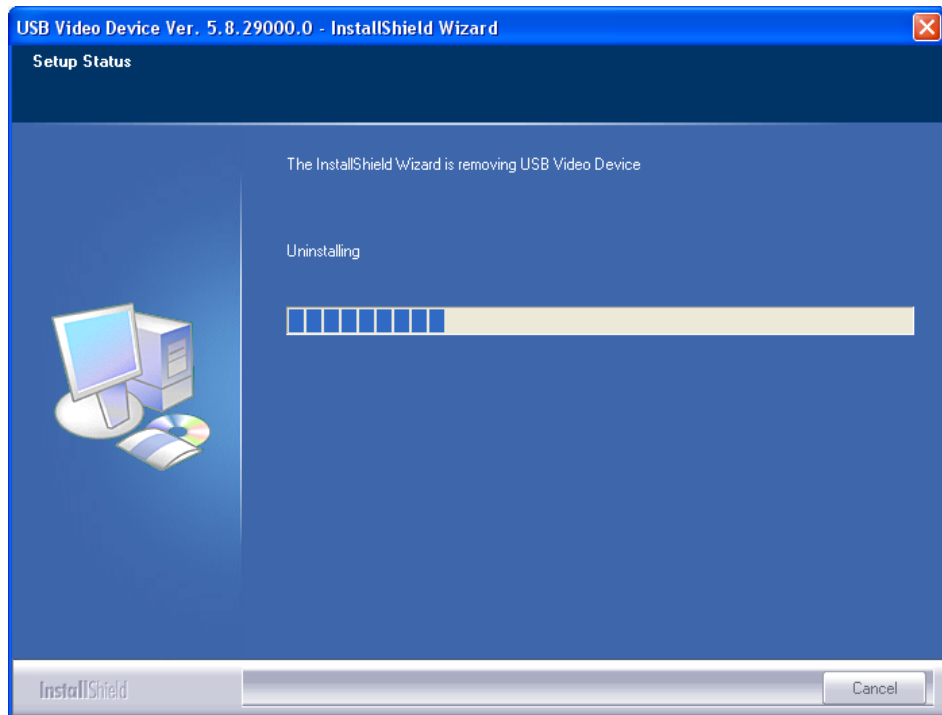


Fig.5-4 in process of uninstall

5.5 Finish uninstall H-CAM USB digital video driver.

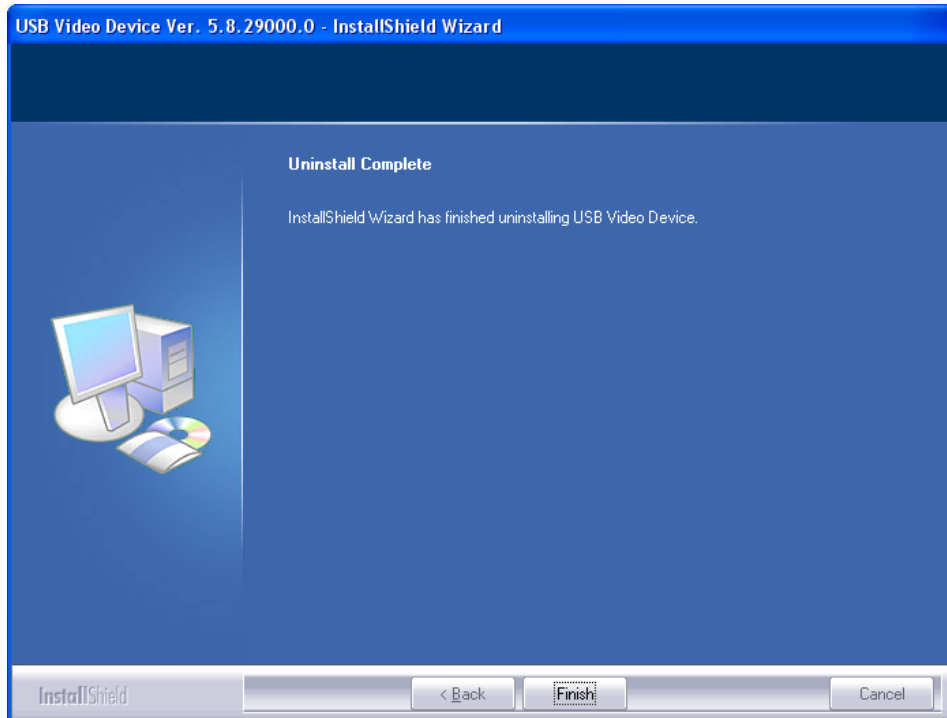


Fig.5-5 Finish uninstall